



EQUIPMENT CALIBRATION RECORD

Equipment Name	UV-Visible Spectrophotometer
Asset Tag / ID	
Type/ Model No.	GENESYS 150
Serial No.	9A5X109012
Location/Department	NARC Analytical Laboratory

Date	Calibrated	Calibration Due Date	Procedure	Remarks	Performed by	Verified by
12-10-2021		12-10-2021	1. Turn on spectrophotometer 2. Select quant in the menu 3. Select if 1 or 2 replicates are needed 4. Selects units in which the sample measurements are displayed 5. Add standards and select curve type 6. Read all the standard series 7. Calibrate button is enabled once concentration values are provided 8. Equation and r-square will be calculated and Measure button is enabled after all standards are measured		ROBERTO T. PISONTE	
add rows if necessary						

Vision:
Mission:

A globally competitive university for science, technology, and environmental conservation.
Development of a highly competitive human resource, cutting-edge scientific knowledge and innovative technologies for sustainable communities and environment.